Metten:

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Measurement:

- 1) Insert the pH-sensor in the liquid.
- 2) Press the **MODE** key to switch-on the instrument and the date **01/01/2020** of the last calibration/reset will be shown in the display followed by the actual pH-value of the liquid that is measured, wait until the value is stable and read of the value.

pH calibration:

Press the MODE key to turn on the instrument

- 1) Insert the pH-sensor in the Nieuwkoop pH-7 calibration liquid.
- 2) Press the **LOG-INC** key for 3 seconds and **ZERO CAL** will appear in the display, alternate by the pH-value.

Note: Wait until the value is stabilised!.

- 3) Press shortly the **SET UP** key and the standard calibration value **6.88** appears in the display (see bottle).
- 4) Press the **SET UP** key for approx. 3 seconds to confirm this value.
- 5) When the value is confirmed **LAST CAL** will appear in the display, followed by **XX / XX / XX** (this is the date of the last calibration), this date can be registered as follows: with the **INC** key you can be change the figure and with **SHIFT** key you can change the position of the cursor, then press the **SET UP** key for 3 seconds to confirm the date and **UPDATE** will appear in the display.

Note: Rinse, clean and dry the sensor before you place the sensor in another liquid.

- 6) Now Insert the pH-sensor in the Nieuwkoop pH-4 calibration liquid.
- 7) Press the **SET UP** key for 3 seconds and **SENS CAL** will appear in the display, alternate by the pH-value.

Note: Wait until the value is stabilised!.

- 8) Press shortly the **SET UP** key and the standard calibration value **4.00** appears in the display (see bottle).
- 9) Press the **SET UP** key for approx. 3 seconds to confirm this value.
- 10) When the value is confirmed **LAST CAL** will appear in the display, followed by **XX / XX / XX** (this is the date of the last calibration), this date can be registered as follows: with the **INC** key you can be change the figure and with **SHIFT** key you can change the position of the cursor, then press the **SET UP** key for 3 seconds to confirm the date and **UPDATE** will appear in the display.

A complete user-manual can be downloaded from the website www.meten.nl

Quick guide Measurement and Calibration pH-meter PH-95

Measurement:

- 1) Insert the pH-sensor in the liquid.
- 2) Press the **MODE** key to switch-on the instrument and the date **01/01/2020** of the last calibration/reset will be shown in the display followed by the actual pH-value of the liquid that is measured, wait until the value is stable and read of the value.

pH calibration:

Press the MODE key to turn on the instrument

- 1) Insert the pH-sensor in the Nieuwkoop pH-7 calibration liquid.
- 2) Press the **LOG-INC** key for 3 seconds and **ZERO CAL** will appear in the display, alternate by the pH-value.

Note: Wait until the value is stabilised!.

- 3) Press shortly the **SET UP** key and the standard calibration value **6.88** appears in the display (see bottle).
- 4) Press the **SET UP** key for approx. 3 seconds to confirm this value.
- 5) When the value is confirmed **LAST CAL** will appear in the display, followed by **XX / XX / XX** (this is the date of the last calibration), this date can be registered as follows: with the **INC** key you can be change the figure and with **SHIFT** key you can change the position of the cursor, then press the **SET UP** key for 3 seconds to confirm the date and **UPDATE** will appear in the display.

Note: Rinse, clean and dry the sensor before you place the sensor in another liquid.

- 6) Now Insert the pH-sensor in the Nieuwkoop pH-4 calibration liquid.
- 7) Press the **SET UP** key for 3 seconds and **SENS CAL** will appear in the display, alternate by the pH-value.

Note: Wait until the value is stabilised!.

- 8) Press shortly the **SET UP** key and the standard calibration value **4.00** appears in the display (see bottle).
- 9) Press the **SET UP** key for approx. 3 seconds to confirm this value.
- 10) When the value is confirmed **LAST CAL** will appear in the display, followed by **XX / XX / XX** (this is the date of the last calibration), this date can be registered as follows: with the **INC** key you can be change the figure and with **SHIFT** key you can change the position of the cursor, then press the **SET UP** key for 3 seconds to confirm the date and **UPDATE** will appear in the display.

A complete user-manual can be downloaded from the website www.meten.nl